



Leica
MICROSYSTEMS

Leica ICC50 E Leica ICC50 W User Manual

Contents

General Notes	
General Notes	4
Symbols Used	5
Safety Instructions	6
Regulatory Compliance Information (Regarding Leica ICC50 W Only)	9
Service	
Care, Maintenance, Contact Persons	12
The Leica ICC50 E or Leica ICC50 W	
Introduction	15
The Basics	16
Leica ICC50 E or Leica ICC50 W: Overview	17
Unpacking	19
Get Ready!	
Preparation for Use	21
Get Set!	
Connection	24
USB Mode – Capturing and Adjusting Images Using a Computer	25
SD Mode – Capturing Images Without a Computer or Without a Wireless Device	27
Adjusting Settings While in SD Mode	29
WiFi Mode – Capturing and Adjusting Images Using a Wireless Device (Only for Leica ICC50 W)	30

Ethernet Mode – Capturing and Adjusting Images Using a Network	31
--	----

The Camera Remote Control (Optional)	
Optional Remote Control	33
Viewing Images and Movies With the Optional Remote Control	34
Calling up the Camera Menu	35
COLOR (Automatic White Balance)	36
COLOR (Manual White Balance)	37
EXPOSURE	38
RESOLUTION	39
SETUP CAMERA (Camera Settings)	40
SETUP USER	41
SETUP ETHERNET	42
SETUP WiFi (Only for Leica ICC50 W)	43
Pairing Cameras With Remotes	44

Go!	
Ready! Set! Go!	46

Specifications	
Specifications	48
Specifications (Only for Leica ICC50 W)	49
Certificates (Only for Leica ICC50 W)	50

General Notes

General Notes

The Leica ICC50 E or Leica ICC50 W cameras include an interactive CD-ROM with all relevant user manuals. Keep it in a safe place, and readily accessible to the user. User manuals and updates are also available for you to download and print from our web site at www.leica-microsystems.com.

Also, look for any available Apps for mobile devices in your iOS or Android App store.

Before installing, operating or using the instruments, read the user manuals listed above. In particular, please observe all safety instructions.

To maintain the unit in its original condition and to ensure safe operation, the user must follow the instructions and warnings contained in these user manuals.

Symbols Used



Read and understand user manual before using this device.



Danger due to hot surface

This symbol warns against touching hot surfaces, e.g. those of light bulbs.

Warning of a danger



This symbol indicates especially important information that must be read and complied with. Failure to comply can cause the following:

- Hazards to personnel
- Functional disturbances or damaged instruments

Warning of hazardous electrical voltage



This symbol indicates especially important information that, if not observed, can cause the following:

- Hazards to personnel
- Functional disturbances or damaged instruments

Important information



This symbol indicates additional information or explanations that intend to provide clarity.

Safety Instructions

Safety concept

Before using your Leica ICC50 E or Leica ICC50 W camera for the first time, please read the "Safety concept" brochure included with your instrument. It contains additional information about handling and care.



Description

The Leica ICC50 E or Leica ICC50 W fulfills the highest requirements for observation and documentation.

Designated use

See "Safety concept" booklet

Non-intended use

See "Safety concept" booklet



The devices and accessories described in this operating manual have been tested for safety and potential hazards. The responsible Leica affiliate must be consulted whenever the device is altered, modified or used in conjunction with non-Leica components that are outside of the scope of this manual.

Unauthorized alterations to the instrument or noncompliant use shall void all rights to any warranty claims!

Location of use

See "Safety concept" booklet



Set up the electrical components at least 10 cm from walls and combustible objects.

Avoid large temperature fluctuations, direct sunlight and vibrations. These conditions can distort measurements and micrographic images.

Requirements for the owner/operator

See "Safety concept" booklet

Ensure that:

The Leica ICC50 E or Leica ICC50 W camera is operated, maintained and repaired by authorized and trained personnel only.

All operators have read, understood and observe this user manual, and particularly the safety instructions.

Safety Instructions (Continued)

Transport

Use the original packaging for shipping or transporting the Leica ICC50 E or Leica ICC50 W camera.

In order to prevent damage from vibrations, disassemble all moving parts that (according to the user manual) can be assembled and disassembled by the customer and pack them separately.

Installation in third-party products

See "Safety concept" booklet

Disposal

See "Safety concept" booklet

Legal regulations

See "Safety concept" booklet

EC Declaration of Conformity

See "Safety concept" booklet

Health risks



Workplaces equipped with microscopes facilitate and improve the viewing task, but they also impose high demands on the eyes and holding muscles of the user. Depending on the duration of uninterrupted work, asthenopia and musculoskeletal problems may occur. For this reason, appropriate measures for reduction of the workload must be taken:

- Optimal arrangement of workplace, work assignments and work flow (changing tasks frequently).
- Thorough training of the personnel, giving consideration to ergonomic and organizational aspects.

Safety Instructions (Continued)

Instrument safety and EMC

Our device has been designed, produced and tested in compliance with

- IEC EN 61010-1: Electrical equipment for measurement, control and laboratory use – Safety – Part one: general requirements
- EN 60950-1: Information technology equipment – Safety – Part one: general requirements
- EN 300328: Electromagnetic compatibility and Radio spectrum Matters (ERM); Wide-band transmission systems; Data transmission equipment operating in the 2,4 GHz ISM band and using wide band modulation techniques; Harmonized EN covering essential requirements under article 3.2 of the R&TTE directive.
- Radio interference suppression in compliance with EN 55011 class B

- Noise immunity in compliance with DIN EN 61326-1.

The device meets the requirements of EU directives

- 2014/35/EU Low voltage directive
- 2014/30/EU EMC directive
- 2011/65/EU RoHS directive

and carries the CE mark.



The device must be disposed of in accordance with 2012/19/EU WEEE directive.

Intended for indoor use only in all EU member states, EFTA states and Switzerland.

Regulatory Compliance Information (Regarding Leica ICC50 W Only)

FCC Requirements for Operation in the United States

FCC radio frequency interference warnings and instructions

This equipment has been tested and found to comply with the limits for a **Class B digital device**, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.

- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

FCC information to user

This product does not contain any user serviceable components and is to be used with approved antennas only.

Any product changes or modifications will invalidate all applicable regulatory certifications and approvals.

FCC guidelines for human exposure

This equipment complies with FCC radiation exposure limits set forth for an uncontrolled environment. This equipment should be installed and operated with minimum distance of 20 cm between the radiator and your body. This transmitter must not be co-located or operating in conjunction with any other antenna or transmitter.

FCC caution

- Any changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate this equipment.
- This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) This device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.
- This device and its antenna(s) must not be co-located or operation in conjunction with any other antenna or transmitter.

Regulatory Compliance Information (Regarding Leica ICC50 W Only) (Continued)

Canada (IC)

This device complies with Industry Canada licence-exempt RSS standard(s). Operation is subject to the following two conditions:

- (1) this device may not cause interference, and
- (2) this device must accept any interference, including interference that may cause undesired operation of the device.

Canada (IC)

Le présent appareil est conforme aux CNR d'Industrie Canada applicables aux appareils radio exempts de licence. L'exploitation est autorisée aux deux conditions suivantes:

- (1) l'appareil ne doit pas produire de brouillage, et
- (2) l'utilisateur de l'appareil doit accepter tout brouillage radioélectrique subi, même si le brouillage est susceptible d'en compromettre le fonctionnement.

Australia and New Zealand Notices

This device equipment complies with the Australian and New Zealand regulatory approvals requirements.



Notice for users in Korea

무선설비의 기기: 제작자 및 설치자는 해당 무선설비가 전파혼신 가능성이 있으므로 인명안전과 관련된 서비스는 할 수 없습니다.

Brasil



Este equipamento opera em caráter secundário, isto é, não tem direito a proteção contra interferência prejudicial, mesmo de estações do mesmo tipo, e não pode causar interferência a sistemas operando em caráter primário.

Service

Care, Maintenance, Contact Persons

General

We hope you enjoy using your high-performance camera. Leica microscope cameras are renowned for their robustness and long service life. Observing the following care and cleaning tips will ensure that even after years and decades, your Leica camera will continue to work as well as it did on the very first day.

Warranty benefits

The guarantee covers all faults in materials and manufacture. It does not, however, cover damage resulting from careless or improper handling.

Care

Keeping all optical components clean is important for maintaining good optical performance.

Cleaning

- Do not use any unsuitable cleaning agents, chemicals or techniques for cleaning.
- Never use chemicals to clean colored surfaces or accessories with rubberized parts. This could damage the surfaces, and specimens could be contaminated by abraded particles.
- Optical surfaces should be cleaned using a lint-free cloth, lens cloth or cotton swab moistened with a commercially available glass cleaner. Do not use alcohol.
- If any optical surface becomes badly coated with dust or dirt brush it off with a camel hair brush before attempting to wipe it clean.

Cleaning polymer components

Some components are made of polymer or are polymer-coated. They are, therefore, pleasant and convenient to handle. The use of unsuitable cleaning agents and techniques can damage polymers.

Care, Maintenance, Contact Persons (Continued)

Protection from dirt

Dust and dirt will affect the quality of your results.

Put the available dust cover over the microscope camera when it will not be used for a long time.

Protect your microscope camera from moisture, fumes and acids and from alkaline, caustic and corrosive materials and keep chemicals away from the instruments.

Protect your microscope camera from oil and grease.

Do not grease guide surfaces or mechanical parts.

Repair, service work

See "Safety concept" booklet

Repairs may only be carried out by Leica Microsystems-trained service technicians.

Only original Leica Microsystems spare parts may be used.

Before opening the instruments, switch off the power and unplug the power cable.

Touching the live circuit can cause injury.

Plugs, optical systems and mechanical parts must not be disassembled or replaced, unless doing so is specifically permitted and described in this manual.

Electrical considerations



Risk of electric shock. Removing the covers of the Leica ICC50 E or Leica ICC50 W camera exposes hazardous voltages, which if contacted, could cause injury or death. Please contact an authorized Leica Microsystems dealer for technical service.

Contact address

If your instrument no longer works perfectly, contact your Leica representative. You can find information on worldwide Leica representatives on the Leica website: www.leica-microsystems.com.

Leica Microsystems (Switzerland) Ltd.
Max Schmidheiny Strasse 201
CH-9435 Heerbrugg (Switzerland)

The Leica ICC50 E or Leica ICC50 W

Introduction

Thank you for purchasing the Leica ICC50 E or Leica ICC50 W by Leica Microsystems. The special design of the Leica ICC50 E or Leica ICC50 W makes it a universal, highly versatile tool for viewing microscopic specimens and capturing still images or video.

Firmware for cameras is constantly being updated. Be sure to keep your camera up-to-date by checking the download section on the camera's web page. There you'll find new firmware and instructions on how to update your camera. You may also contact your Leica representative for help.

The Basics

i Your Leica HD camera has many possibilities for various applications and conditions of use. Below are some basic concepts to help you get the most performance for your application.

Power

Leica ICC50 E or Leica ICC50 W – 5 V/2 A

Power can be provided to the camera in the following ways:

- USB power connector on the microscope stand
- External USB power supply
- USB connectors PC

Modes of use

1. SD mode

- capture onto SD card
- use remote control

2. WiFi mode (only available for ICC50 W)

- Wireless connection to wireless device (tablet, phone, PC)

3. USB mode


- Cabled connection to PC

4. Ethernet mode

- Cabled connection to external network

5. Capture button

- Captures image onto the SD card

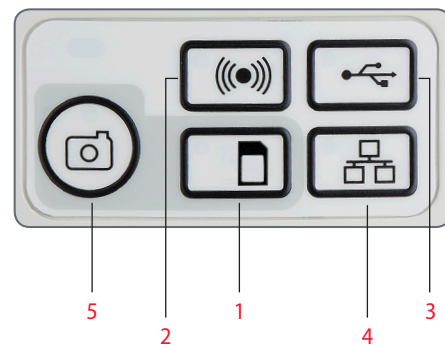

i Note direct connection to an HD display is possible in all modes. Aspect ratio is 16:9 except in USB mode where the aspect ratio is 4:3.

Leica ICC50 E or Leica ICC50 W: Overview



1. LED status indicator

- *Blinking red means starting up*
- *Green means ready*

2. Capture button for SD card capture

3. Mode selection panel


4. SD card slot

5. Power on/off button


6. Service button



Leica ICC50 E or Leica ICC50 W: Overview (Continued)

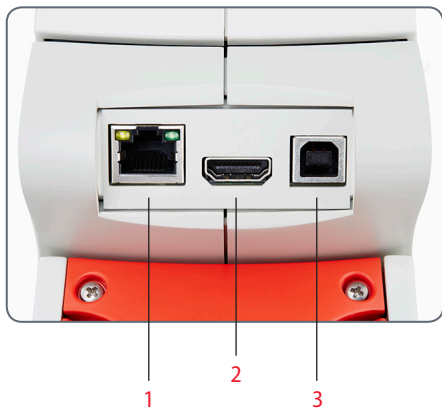


1. *Ethernet socket for connecting an ethernet cable (cable not provided)*
2. *HDMI for connecting the provided HDMI cable*
3. *USB for connecting the provided USB cable*

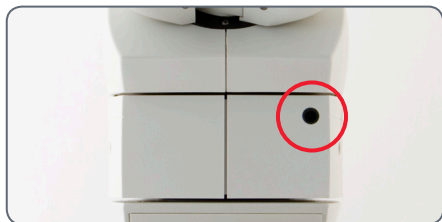

Front

Infrared window: Receives the signal from the optional remote control.

Unpacking

 Please do not discard any of the packing materials. They should be used for safely storing and transporting the instrument should the need arise.



The contents include the camera module, CD with imaging software, HDMI cable, and USB cable.



If you ordered the Optional USB power supply you have also received a power supply.




If you ordered the handheld remote, you received the remote with batteries.


Get Ready!


Preparation for Use

Tools used

- Allen key (provided with microscope)

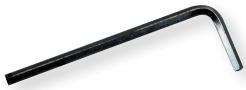

 Do not insert anything except the SD memory card into the card slot – danger of short circuit and electric shock!

 The Leica ICC50 E or Leica ICC50 W camera module fits between the microscope stand and the microscope's viewing tube.

Preparation for use

1. If the microscope's viewing tube is in place, remove it.

Attach the Leica camera to the microscope stand:

2. Slightly unscrew the set screw (located on top of the stand) with the allen key tool provided with the microscope.



3. Insert the dovetail (see the underside of the Leica ICC50 E or Leica ICC50 W) into the stand support.




4. Tighten the set screw firmly without forcing it.



Preparation for Use (Continued)


5. Attach the microscope viewing tube onto the Leica ICC50 E or Leica ICC50 W using the set screw, which is located on the left side of the camera module.

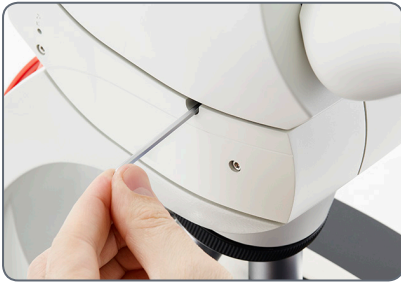

Get Set!

Connection

Connect the camera to a power source


 Note the Leica ICC50 E or Leica ICC50 W requires 5 V/2 A for power, therefore, a proper power supply must be used.

Connect to a power source by one of the following methods:

1. Via provided USB cable to the USB connector on back of the microscope stand.



2. Via provided USB cable to external USB power supply.

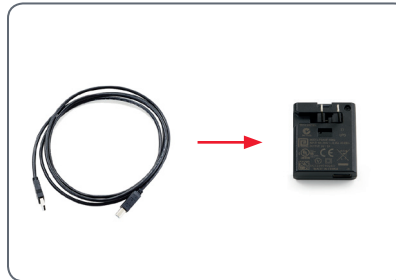

3. Via provided USB cable to USB connector of a PC. Camera should be in USB mode.




USB Mode – Capturing and Adjusting Images Using a Computer

i 10× eyepieces have a fixed circular field of view of 18 mm or more diameter. The Leica ICC50 E or Leica ICC50 W's field of view, however, is rectangular with a diagonal of 16.4 mm to ensure an even field on the digital image. As a result, the image field of the camera is smaller than the object field seen through the eyepieces.

Installing and connecting

1. Insert the provided CD into your computer, and follow the instructions to load the software.

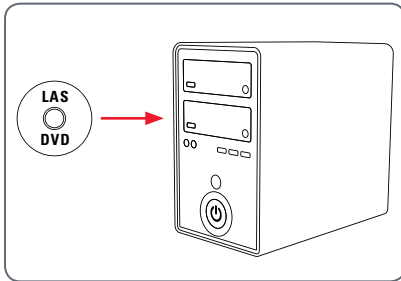

2. Insert the provided USB cable into the camera's USB port, and attach the other end of the cable to a USB 2.0 port of your computer.



3. Press the red power button on the camera.



4. The status light will blink red. Wait until the status light is green and one of the mode button's is also green. This may take approximately 30 seconds.





USB Mode – Capturing and Adjusting Images Using a Computer (Continued)


5. If a mode button other than USB mode is lit green, then press the USB mode button and wait until both the status LED and the USB mode button is green.



6. You are ready to launch the Leica software. Follow the software instructions to adjust and capture an image.
7. While in USB mode, you can connect the provided HDMI cable from the camera to an HD monitor or HD projector to share the image.

 Note: You won't be able to write or read the SD card, if the camera is connected via USB cable with your PC.

 Note the image aspect ratio while in USB mode is 4:3. To avoid distortion if your projector or monitor is set to a 16:9 aspect ratio, you will need to adjust the aspect ratio of the projector or monitor to 4:3 using the controls of the projector or monitor.

 Note the image resolution on the HD projector or monitor is limited to the live image resolution setting in the software.




4:3 aspect ratio



4:3 aspect ratio image on a screen set to an aspect ratio of 16:9

SD Mode – Capturing Images Without a Computer or Without a Wireless Device

 In order to use this mode, you must have an SD card (not provided) with free memory on it. If the SD memory card is full, the capture button LED will not light green and you will not be able to capture any more images.

Attaching the power cable

1. Be sure the camera is connected to a power source – see power source section on [page 24](#).
2. Press the red power button.



3. The status light will blink red. Wait until the status light is green and one of the mode button's is also green. This may take approximately 30 seconds.



4. If a mode button other than SD mode is lit green, then press the SD mode button and wait until both the status LED and the SD mode button is green.



SD Mode – Capturing Images Without a Computer or Without a Wireless Device (Continued)

5. Insert the SD memory card into the slot on the side of the Leica ICC50 E or Leica ICC50 W until the card clicks into place.



The capture button LED will turn green. The Leica ICC50 E or Leica ICC50 W is now ready to capture images onto the SD memory card.



Capturing an image



It is critical to properly set the microscope eyepieces' diopters to ensure that the images on the SD memory card are in focus when the image is in focus through the microscope's eyepieces. Refer to the microscope's user manual for proper diopter setting. Focus the microscope at high magnification, then change to the magnification at which you wish to capture the image. Do not refocus the microscope. You can also use the HDMI output and an HD display (not provided) as a focusing aid.


6. To capture an image onto the SD memory card, lightly press the capture button, located on the side of the Leica ICC50 E or Leica ICC50 W camera.



- You will hear a beep to acknowledge the button was pressed.
- The status LED flashes red while the image is being captured.
- You will hear another beep when the Leica ICC50 E or Leica ICC50 W has completed capturing the image and the status LED will return to steady green.


The image is now saved on the SD card. Images which are on the SD card can be viewed using the optional remote control. See "Camera Remote Control" on [page 32](#).

7. While in SD mode, you can connect the provided HDMI cable from the camera to an HD monitor or HD projector to share the image.

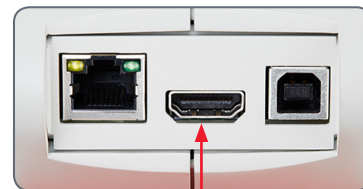

Adjusting Settings While in SD Mode

Setting a new white balance

 Make sure that there is no sample in the beam path and that a standard illumination is available when you reset the camera.


Press and hold the capture button for 5 seconds. Wait for the beep. A new white balance is set.



Return to factory settings

Press and hold the service button for 5 seconds. Wait for the beep. The camera will now reset to factory settings.



 It is recommended to set a new white balance each time you change the type or color temperature of the illumination. This is especially needed when you work with halogen light which can be turned from low intensity yellow to a high intensity blue light. The new white balance setting will generate a neutral background for both illuminations.

WiFi Mode – Capturing and Adjusting Images Using a Wireless Device (Only for Leica ICC50 W)


1. Be sure the camera is connected to a power source – see power source section on [page 24](#).



2. Press the red power button.
3. The status light will blink red. Wait until the status light is green and one of the mode button's is also green. This may take approximately 30 seconds.




4. If a mode button other than WiFi mode is lit green, then press the WiFi button and wait until both the status LED and the WiFi mode button is green.

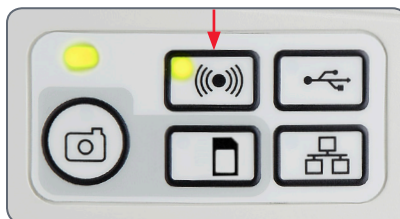

5. Follow the directions of your WiFi viewing device to connect to the internal WiFi network of the camera.
WiFi Password: leicamicro
6. Open the appropriate app on your mobile device or appropriate software on your wireless enabled PC for viewing, capturing, and adjusting the camera image.

Ethernet Mode – Capturing and Adjusting Images Using a Network

1. Connect one end of an ethernet cable (not provided) to the camera and the other end of the cable to the ethernet connection in your facility.

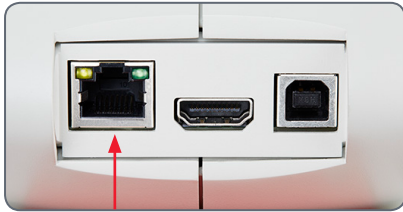

2. Be sure the camera is connected to a power source – see power source section on [page 24](#).
3. Press the red power button.



4. The status light will blink red. Wait until the status light is green and one of the mode button's is also green. This may take approximately 30 seconds.



5. If a mode button other than ethernet mode is lit green, then press the ethernet button and wait until both the status LED and the ethernet LED mode button is green.



6. Follow the directions of your network compatible device to connect to your facilities network and select the camera.
7. Open the appropriate app on your mobile device or appropriate software on your wireless enabled PC for viewing, capturing, and adjusting the camera image.

The Camera Remote Control (Optional)



The remote control is designed to work only when the camera is in SD mode and when the camera is used in combination with an HD display.

The only function which works in any mode is the "Info" function which shows details about the camera.

Optional Remote Control




The remote control can only work in SD mode except for the "i" button which will display information about the camera on an HD display. The remote must be pointed at the IR receiver of the camera without obstructions.







1. Start/stop video recording
The max. recording time of a movie is 23.00 min. After this time you have to restart the recording again.
2. Save still image to SD card
3. Pause/play movie clip
4. View files on SD card
5. Move to previous image on SD card
6. Move to next image on SD card
7. Brightness adjustment up or down
8. Pairing button
9. Show/hide information box
10. White balance set button
11. Call up the camera menu
12. Arrow keys for navigation
13. OK to confirm

Viewing Images and Movies With the Optional Remote Control

 If you have purchased the optional remote, the Leica HD camera shows images and movie clips directly on an HD monitor.


1. Thumbnail view  (first press)

- up/down/left/right ► select thumbnail
-  Show image in full screen or playback movie clip
- next image  / previous  image
-  ► Play movie clip
- left/right ► fast forward / rewind

2. To return to live view press .



Playback / fullscreen view

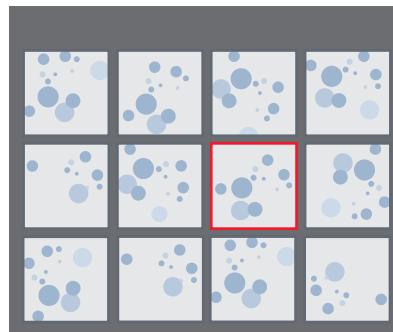








Thumbnail view

Calling up the Camera Menu

 The remote control can only work when the camera is in SD mode.




Opening and closing the camera menu


1. Point the remote control towards the camera.
2. Press the  key on the remote control to show the camera menus on the monitor.
3. Press the , ,  and  buttons to select a menu item.
4. Press the  button to confirm a menu item.
5. Press the  button again to hide the camera menus.



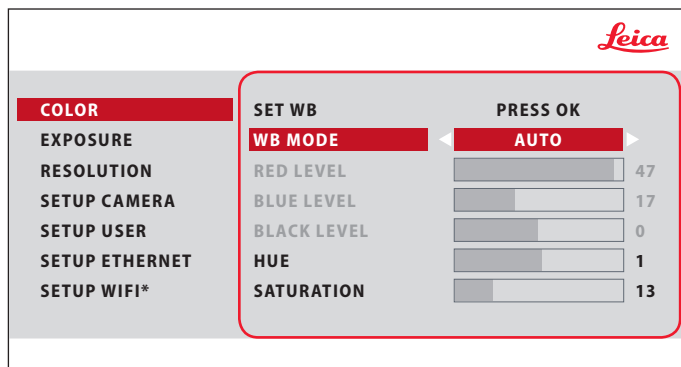

COLOR (Automatic White Balance)




* SETUP WIFI only for ICC50 W


 The "COLOR" function makes it possible to adapt the camera chip to the ambient light so that color-neutral images can be acquired.

Enabling automatic white balance

1. Press the  button on the remote control.
2. Call up the "COLOR" entry.
3. Set the value for "WB MODE" to "AUTO".
4. Press the  button to exit the menu.

 The values for "RED" and "BLUE" cannot be adjusted if the "WB MODE" setting is set to "AUTO".

COLOR (Manual White Balance)

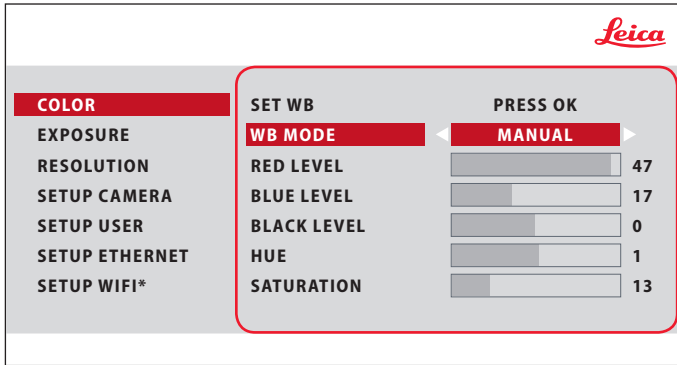



* SETUP WIFI only for ICC50 W




The "COLOR" function makes it possible to adapt the camera chip to the ambient light so that color-neutral images can be acquired.


Adjusting the white balance manually (recommended)


1. Move the specimen out of the field of view so that only the illumination is seen.
2. Press the  button on the remote control.
3. Call up the "COLOR" entry.
4. Set the "WB MODE" to "MANUAL".
5. If necessary, correct the values for "RED" and "BLUE" until the image shows a uniform gray area.
6. Press the  button to exit the menu.

EXPOSURE

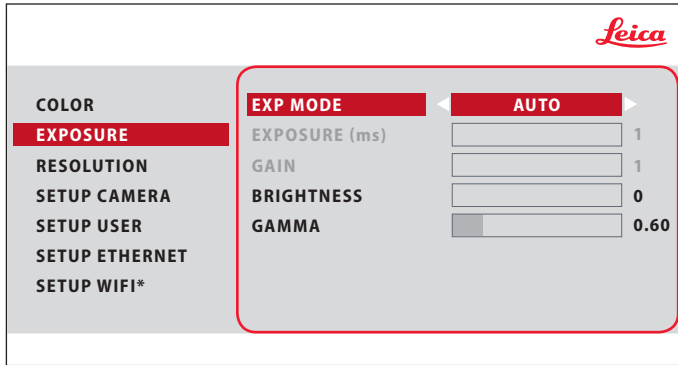


Automatic exposure

1. Press the  button on the remote control.
2. Call up the command for automatic exposure (see diagram at the left).
3. Set the value for "EXP MODE" to "AUTO".
4. Correct the values for "BRIGHTNESS" and "GAMMA" until you obtain the desired results for the image.

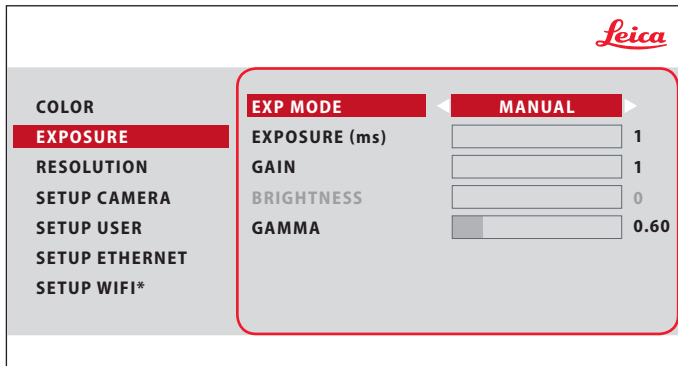


Manual exposure


1. Press the  button on the remote control.
2. Call up the command for manual exposure (see diagram at the left).
3. Set the value for "EXP MODE" to "MANUAL".
4. Correct the values for "EXPOSURE", "GAIN" and "GAMMA" until you obtain the desired results for the image.

* SETUP WIFI only for ICC50 W

RESOLUTION

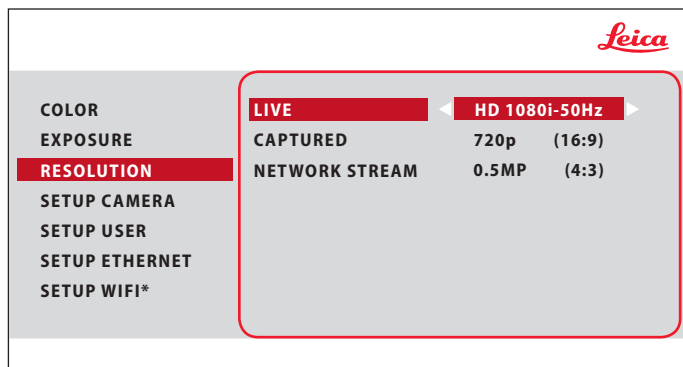

* SETUP WIFI only for ICC50 W

 The "RESOLUTION" menu enables you to define the resolution for the live image, still image capture and network stream individually. This provides optimum results in every situation. Video capture on SD card will always be 1,920 × 1,080 resolution.

LIVE (resolution of the live image)

Defines the resolution of the live image. After changing the live resolution, you will see a message on the HD screen that you can accept the new resolution by pressing "OK" otherwise, the live resolution will return to the previous setting.


CAPTURED (resolution of the captured image)

Defines the resolution of the captured image on the SD card.


NETWORK STREAM

Indicates the resolution of the live image when in WiFi or Ethernet mode.

SETUP CAMERA (Camera Settings)

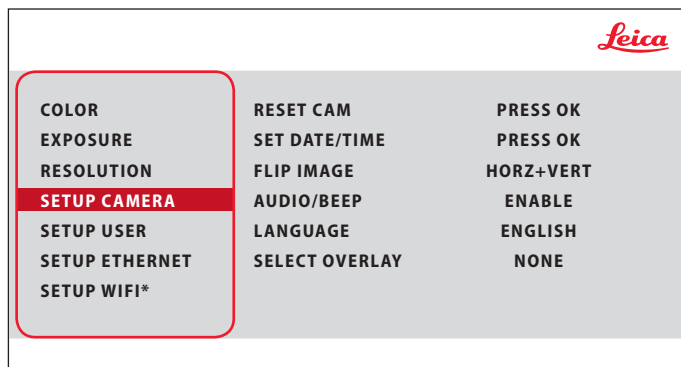

* SETUP WIFI only for ICC50 W

 The camera settings enable you to control internal parameters such as the date or time, reset the camera, etc.

RESET CAMERA

This command resets the camera to the factory settings. All user information (such as white balance, resolution etc.) is reset.

SET DATE TIME

This command sets the date and time as well as the format for displaying them. The following formats are available for selection: YYYYMMDD – DDMMYY – MMDDYY

FLIP IMAGE

Select to flip image 180°.


AUDIO/BEEP


You can decide to have a tone from the camera for some functions.

LANGUAGE


Select the language of the On Screen Display.

SELECTING IMAGE OVERLAYS (SELECT OVERLAY)

Press ◀ ▶ to select another overlay. Press  to activate the selected overlay and leave the menu.

 If you want to quickly display other overlays, press when the live image is displayed ▼ ▲ to show the next or previous respective overlay.

SETUP USER


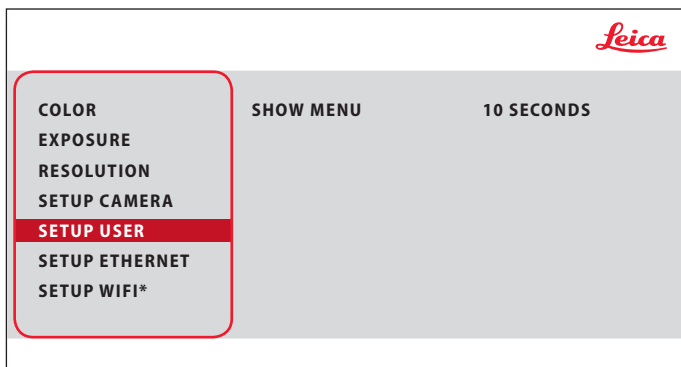


SETUP USER


Here you can decide to have the menu remain on screen for 10 seconds or until you press the menu button again.

* SETUP WIFI only for ICC50 W

SETUP ETHERNET

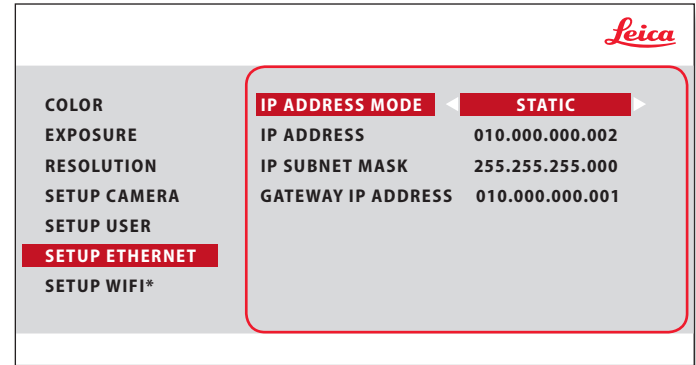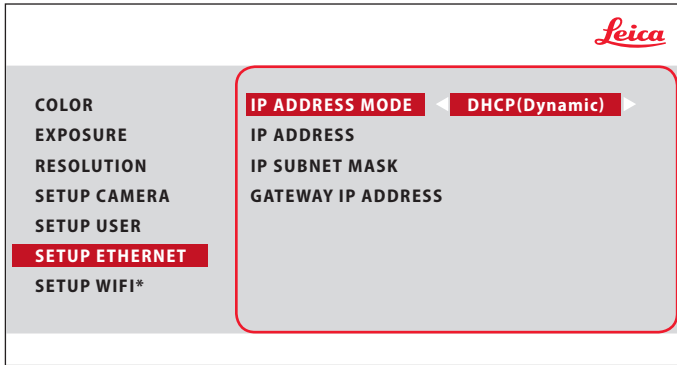

* SETUP WIFI only for ICC50 W

 This menu allows you to customize parameters relevant to working in Ethernet mode. Consult with your local IT expert for information relating to your local network.


IP ADDRESS

You can decide if you want the IP address to be set by your network (Dynamic) or by you using the onscreen keyboard (Static).


If you selected "STATIC" for the IP address you can input the following using the onscreen keyboard:

- IP address
- IP subnet mask
- Gateway IP address

SETUP WiFi (Only for Leica ICC50 W)

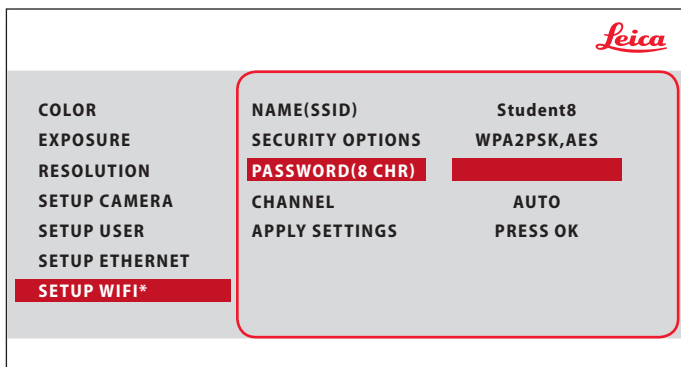

* SETUP WIFI only for ICC50 W

 This menu allows you to customize parameters relevant to working in WiFi mode.

NAME (SSID)

This is the camera name which will appear. You can change the name using the onscreen keyboard.

SECURITY OPTIONS – NONE

This means there is no password.

SECURITY OPTIONS – WPA2-PSK AES

This allows you to create a password.

PASSWORD

Create a password using the onscreen keyboard. Default password is "leicamicro".

CHANNEL


This allows the WiFi channel to be set automatically or by the user.


APPLY SETTINGS


After making any changes to the WiFi settings, you must go to "APPLY SETTINGS" and press "OK" on the remote to implement the changes.


Pairing Cameras With Remotes


Pairing

 The Leica ICC50 E or Leica ICC50 W and the remote control can be paired and then only respond to each other. This can be helpful when using multiple cameras and remote controls.

1. Press the  button to start or end the process.

 In order to achieve successful pairing and avoid pairing by mistake, the second step must be performed within 4 seconds.

 If a "timeout" is displayed after 4 seconds, press the "Pair" button again to start the process.

2. Press and hold the button on your remote you want to define as the pairing button until a corresponding confirmation is displayed on the HD monitor. All of the buttons except the  button can be used for this.

3. The microscope only responds to the remote control commands from this remote control.

Resetting to factory setting

1. Press the button  to start the process.
2. Press the button  until a corresponding confirmation is displayed on the HD monitor.

Go!

Ready! Set! Go!

Now all you need to do is properly set up your microscope, capture your images, and enjoy!

Specifications

Due to the technology of the camera, the specifications are changing regularly. Please visit our website www.leica-microsystems.com/education for the latest brochure downloads to find the up to date specifications.

Specifications

Operating environment

Feature	Specification
Temperature	
Operating	+10°C to +40°C (50°F – 104°F)
Storage	-20°C to +55°C (-4°F – 131°F)
Relative humidity (maximum)	
Operating/Storage	10 % to 90 % (noncondensing)
Altitude (maximum)	
Operating	0 m to 2,000 m (0 ft to 6,561 ft)
Nonoperating	0 m to 10,668 m (0 ft to 35,000 ft)
Pollution Degree	
Rating	2
Weight	0.7 kg




Ports and connectors

Feature	Specification
USB	one standard USB 2.0 (Typ B)
Video	one standard HDMI connector (Typ A)
Network adapter	one RJ-45 connector
Memory card reader	supports up to 32 GB
Power supply	via USB cable to external power supply, microscope stand or PC use USB 2.0 or higher (do not use lower USB version for power input)
Power requirement	max. 10 W

Specifications (Only for Leica ICC50 W)

Feature	Specification
Radio technology	IEEE 802.11n (HT20)
Operating frequency	2,412 – 2,472 MHz (20 MHz)
Transmitted power	26.3 mW /14.2 dBm (20 MHz) e.i.r.p

Certificates (Only for Leica ICC50 W)

Declaration of Conformity Konformitätserklärung Déclaration de Conformité		
We / Wir / Nous		
Leica Microsystems (Schweiz) AG Industry Division Max Schmidheiny-Strasse 201 CH-9435 Heerbrugg		
declare under our sole responsibility that the product erklären in alleiniger Verantwortung, dass das Produkt déclare sous notre seule responsabilité que le produit		
13 613 735 ICC50 W Kamera Modul		
To which this declaration relates is in conformity with the following standards Auf das sich diese Erklärung bezieht, mit den folgenden Normen übereinstimmt Auquel se réfère cette déclaration est conforme aux normes		
EN 61010-1:2010 EN60950-1:2006+A11:2009+A11:2010+A12:2011+A2:2013 EN 61326-1:2013 EN 55011:2016 EN62311:2008 EN301489-17 V3.1.1:2017 EN301489-1 V2.1.1:2017 EN300328 V2.1.1:2016 EN 50581:2012		
Following the provisions of directive(s) gemäss den Bestimmungen der Richtlinie(n) conformément aux dispositions de(s) directive(s)		
Radio Equipment Directive (RED) – 2014/53/EU Restriction of the use of certain hazardous substances in electrical and electronic equipment (RoHS) 2011/65/EU		
Place, date, Ort, Datum/ lieu, date	Name and function/ Name und Funktion/ nom et fonction	
Heerbrugg, July 14th, 2017		
	Michael Stroehle Leica Microsystems (Schweiz) AG Global Manager Regulatory/Quality engineering	



Leica Microsystems (Schweiz) AG · Max-Schmidheiny-Strasse 201 · 9435 Heerbrugg, Switzerland
T +41 71 726 34 34 · F +41 71 726 34 44

www.leica-microsystems.com

CONNECT
WITH US!

